

# OneStep PAGE gel

RUO Research Use Only

REF 16171



## INTRODUCTION

OneStep PAGE gel enables researchers to cast, polymerize and run a mini-gel in less than 15 minutes. Provided as a novel, ready-to-pour SDS polyacrylamide solution, it polymerizes into a unique support matrix for the electrophoretic separation of denatured proteins. The proprietary chemistry of OneStep PAGE gel eliminates the need for a stacking gel while providing sharp band resolution over a wide molecular weight range. The gels are fully compatible with all standard electrophoresis equipment, SDS-PAGE staining procedures and downstream application including 2D electrophoresis, Western blot transfer, protein sequencing and MALDI analysis.

## KIT CONTENTS

Label	Contain
OneStep PAGE Solution	125 ml
OneStep PAGE 2x Buffer	125 ml

## STORAGE AND STABILITY

- Storage condition: The product remains active for at least 2 year stored at 4°C in an unopened container.

## FEATURE

- Use general loading buffer and running buffer (Tris-Glycine-SDS buffer)
- Faster gel casting with no stacking gel required
- Broad range of separation
- High resolution of protein bands

Concentration	Protein size
7.5%	20 – 300 kDa
10%	10 – 200 kDa
12%	3.5 – 100 kDa
15%	2.5 – 100 kDa

## APPLICATIONS

- SDS-PAGE electrophoresis

## PROTOCOL

- Assemble clean glass plate and spacer.
- Mix OneStep PAGE solution, 2X buffer, 10% APS and TEMED in the tube.
- Pour the mixed solution into gel casting.
- Insert comb and wait 10~20min for polymerization.
- Remove comb and load sample.
- Run the gel 120V – 150V constant voltage for 60~90min or your protocol.

Contents	SDS-PAGE Gel Percentage (%)			
	7.5%	10%	12%	15%
D.W	2.5ml	1.67ml	1ml	-
OneStep PAGE Solution	2.5ml	3.3ml	4ml	5ml
OneStep PAGE 2x Buffer	5ml	5ml	5ml	5ml
10% Ammonium persulfate	60ul	60ul	60ul	60ul
TEMED	6ul	6ul	6ul	6ul

## ORDERING INFORMATION

Product Name	Amount	Cat. No.
WEST-ZOL™ plus Western Blot Detection System	200ml	16024
PRO-PREP™ Protein Extraction Solution (C/T)	100ml	17081
PRO-MEASURE™ Protein Measurement Solution	100ml	21011
Gang-Nam-STAIN™ Prestained Protein Ladder	250ul	24052
Ponceau S Solution	1L	31192
SenSafe™ Fast Protein Staining Solution	1L	31194

Technical support : +82-505-550-5600

Review date : 2015. 06. 30